

YTzol Pure RNA

Cat No:YT9066	Size: 20 ml
YT9065	Size: 50ml
YT9064	Size:100 ml

Store at 2-8°C, protect from light. YTzol reagent is stable for up to 12 months, when properly stored.

Reagents required, but not supplied:

- Chloroform
- Isopropyl alcohol
- 75% Ethanol (in RNA-free water)
- RNase-free water or 0.5% SDS solution [To prepare RNase-free water, draw water into RNase-free glass bottles. Add diethylpyrocarbonate (DEPC) to 0.01% (v/v). Let stand overnight and autoclave. The SDS solution must be prepared using DEPC-treated, autoclaved water.]

WARNING: Toxic in contact with skin and if swallowed. Causes burns. After contact with skin, wash immediately with plenty of water. If you feel unwell, seek medical advice (show label where possible). Phenol (108-95-2) and other Components (NJTSRN 80100437-5000p).

Precautions for Preventing RNase Contamination

RNases can be introduced accidentally into the RNA preparation at any point in the isolation procedure through improper technique. Because RNase activity is difficult to inhibit, it is essential to prevent its introduction. The following guidelines should be observed when working with RNA.

- Always wear disposable gloves. Skin often contains bacteria and molds that can contaminate an RNA preparation and be a source of RNases. Practice good microbiological technique to prevent microbial contamination.
- Use sterile, disposable plasticware and automatic pipettes reserved for RNA work to prevent cross-contamination with RNases from shared equipment. For example, a laboratory that is using RNA probes will likely be using RNase A or T1 to reduce background on filters, and any nondisposable items (such as automatic pipettes) can be rich sources of RNases.
- In the presence of YTzol Reagent, RNA is protected from RNase contamination. Downstream sample handling requires that nondisposable glassware or plasticware be RNase-free. Glass items can be baked at 150°C for 4 hours, and plastic items can be soaked for 10 minutes in 0.5 M NaOH, rinsed thoroughly with water, and autoclaved.

Recommended volume of YTzol on different starting materials :

10 cm ² adherent cells	1 ml
10 ⁷ suspension cells	1-2 ml
100 ul white cells	2 ml
50-100 mg ordinary tissue	1 ml
50-100 mg special tissue(live, spleen, bone or cartilage)	2 ml
15-100 mg plant tissue	1 ml

Protocol

1. HOMOGENIZATION:

Use YTzol Reagent to prepare lysates from various sample types as described below.

a. Tissues: from animal or plant (either fresh or frozen at -70°C until use) can be processed by freezing with liquid nitrogen and grinding into a fine powder using a mortar and pestle. Homogenize tissue samples in 1 ml YTzol Reagent per 50–100 mg tissue using a tissue homogenizer or rotor-stator.

b. Adherent Cells: Lyse cells directly in a culture dish by adding 1 ml of YTzol Reagent to a 3.5 cm diameter dish, and passing the cell lysate several times through a pipette. The amount of YTzol Reagent added is based on the area of the culture dish (1 ml per 10 cm^2) and not on the number of cells present. An insufficient amount of YTzol Reagent may result in contamination of the isolated RNA with DNA.

c. Suspension Cells: Pellet cells by centrifugation. Lyse cells in YTzol Reagent by repetitive pipetting. Use 1 ml of the reagent per $5\text{-}10 \times 10^6$ of animal, plant or yeast cells, or per 1×10^7 bacterial cells. Washing cells before addition of YTzol Reagent should be avoided as this increases the possibility of mRNA degradation. Disruption of some yeast and bacterial cells may require the use of a homogenizer. **OPTIONAL:** An additional isolation step may be required for samples with high content of proteins, fat, polysaccharides or extracellular material such as muscles, fat tissue, and tuberous parts of plants. Following homogenization, remove insoluble material from the homogenate by centrifugation at $12,000 \times g$ for 10 minutes at 2 to 8°C . The resulting pellet contains extracellular membranes, polysaccharides, and high molecular weight DNA, while the supernatant contains RNA. In samples from fat tissue, an excess of fat collects as a top layer which should be removed. In each case, transfer the cleared homogenate solution to a fresh tube and proceed with chloroform addition and phase separation as described.

2. PHASE SEPARATION :

Incubate the homogenized samples for 5 minutes at 15 to 30°C to permit the complete dissociation of nucleoprotein complexes. Add 0.2 ml of chloroform per 1 ml of YTzol Reagent. Cap sample tubes securely. Shake tubes vigorously by hand for 15 seconds and incubate them at room temperature for 2 to 3 minutes. Centrifuge the samples at $12,000 \times g$ for 5-10 minutes at 4°C . Following centrifugation, the mixture separates into a lower yellow, phenol-chloroform phase, an interphase, and a colorless upper aqueous phase. RNA remains exclusively in the aqueous phase. The volume of the aqueous phase is about 60% of the volume of YTzol Reagent used for homogenization.

3. RNA PRECIPITATION

Transfer the aqueous phase to a fresh tube, and save the organic phase if isolation of DNA or protein is desired. Precipitate the RNA from the aqueous phase by mixing with isopropyl alcohol. Use 0.5 ml of isopropyl alcohol per 1 ml of YTzol Reagent used for the initial homogenization. Mix well by inverting the tube several times. Incubate samples at 15 to 30°C for 10 minutes and centrifuge at $12,000 \times g$ for 5-10 minutes at 4°C . The RNA precipitate, often invisible before centrifugation, forms a gel-like pellet on the side and bottom of the tube.

4. RNA WASH

Discard the supernatant, add 1 ml 75% Ethanol (in RNA-free water), do not stir the precipitate, gently inverting the tube several times to wash the tube, centrifuge at $12,000 \times g$ for 2 min at 4°C , discard the ethanol, repeat the step again.

5. REDISSOLVING THE RNA

At the end of the procedure, briefly dry the RNA pellet (air-dry or vacuum-dry for 5-10 minutes). Do not dry the RNA by centrifugation under vacuum. It is important not to let the RNA pellet dry completely as this will greatly decrease its solubility. Partially dissolved RNA samples have an A_{260}/A_{280} ratio < 1.6 . Dissolve RNA in RNase-free water or 0.5% SDS solution by passing the solution a few times through a pipette tip, and incubating for 10 minutes at 55 to 60°C . (Avoid SDS when RNA will be used in subsequent enzymatic reactions.) RNA can also be redissolved in 100% formamide (deionized) and stored at -70°C .